

Data sheet

This data sheet provided for information only

DiaTaq^{RED} DNA POLYMERASE

RESEARCH USE ONLY

Recombinant DiaTagRED DNA Polymerase

(Deoxynucleosidetriphosphate: DNA Deoxynucleosidyltransferase E.C. 2.7.7.7.)

SOURCE:

Thermus Augaticus YT1 strain

Cat.No.	Pack	Conc.
DTR-250	250 Units	5 U/μl
DTR-1000	1000 Units	5 U/μl

Stability:

Shelf life 24 months if store at -20°C

UNIT DEFINITION

One unit defined as the amount of enzyme that incorporates 10 nmoles of dNTP's into acid-insoluble form in 30 minutes at 74°C under assay conditions.

DESCRIPTION

DiaTaqRed DNA Polymerase is a thermostable 94 kDa DNA Polymerase purified from E.coli PVG-AI recombinant strain expressing Thermus aquatiqus polymerase gene.

The enzyme catalyzes polymerization of nucleotides into duplex DNA in the 5'-3' direction in presence of Mg²⁺ ions. The enzyme is stained in red. The stain facilitates the visual control of mixing the enzyme with other components in reaction. The stain doesn't influence on enzymatic activity of polymerase and other general characteristics

STORAGE AND DILUTION BUFFER:

20mM Tris-HCL (pH 8.0);100mM KCL;0.1mM EDTA; 1mM DTT; 50 glycerol; 0.5% Nonidet P-40; 0.5% Tween-20

AMPLIFICATION BUFFERS:

10X NH4-buffer: 166mM (NH₄)₂SO₄; 670mM Tris-HCL (pH 8.8 at 25°C); 0.1% Tween-20.

"Enhanced" NH4-buffer: 166mM (NH₄)₂SO₄; 670mM Tris-HCL (pH 8.8 at 25°C); 0.1% Tween-20; stabilizers/enhancer

ASSOCIATED ACTIVITIES:

Endonuclease and exonuclease activities were not detectible after 2 and 1 hours incubation, respectively, of 1 µg lambda DNA and 0.22 µg of EcoR I digested lambda DNA, respectively, at $72^{\bar{0}}$ C in the presence of 15-20 units of **DiaTag**^{Red} DNA polymerase

APPLICATIONS:

- Routine PCR
- Primer extension
- RT-PCR

STORAGE CONDITIONS:

Store **DiaTag^{Red}** DNA Polymerase at -20^oC for long term storage

SHIPPING CONDITIONS:

Should be shipped at ambient temperature. For long distance shipments preferably in **Blue Ice**

DIALAT Ltd., 117 149, Moscow, Russia, Sympheropolsky blv.8, Tel. +499 613 2052, Fax. +499 613 4818 e-mail: info@dialat.ru web: www.dialat.ru